



Diamond Stick





Dose the stick in a spiral formation while your disk is running **1**

Alternatively you can dose the diamond stick in a star formation on the disk before running **2**

Dose lubricant $\frac{1}{3}$ outwards on the cloth **3**

Lubricate often with small amounts **4**

You can quickstart a new cloth by using the Diamond Stick.

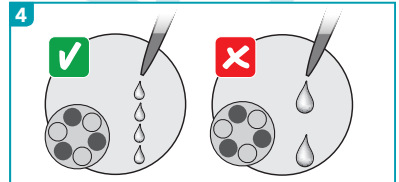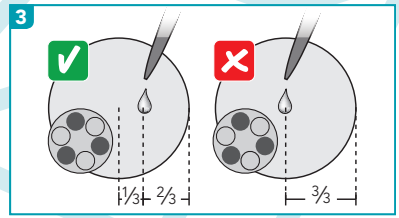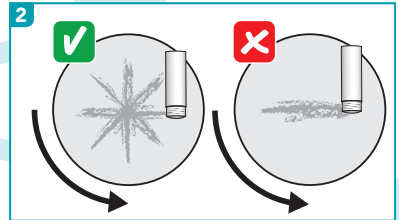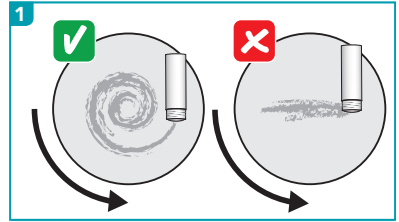

Mono Crystalline
9 μ m - 1 μ m
20 ml



Poly Crystalline
9 μ m - 0.25 μ m
20 ml

Diamantstick





Dosieren Sie den Stick in Spiralform, während die Scheibe läuft **1**

Alternativ können Sie bevor die Scheibe läuft den Stick in Sternform dosieren **2**

Das Schmiermittel sollte im äußeren Drittel der Scheibe aufgetragen werden **3**

Verwenden Sie das Schmiermittel öfters jedoch in geringen Mengen **4**

Mit dem Diamantstick ist das Poliertuch sofort Einsatzbereit.

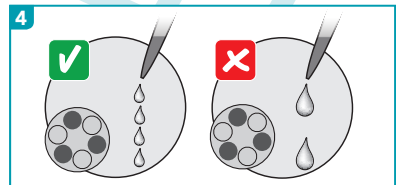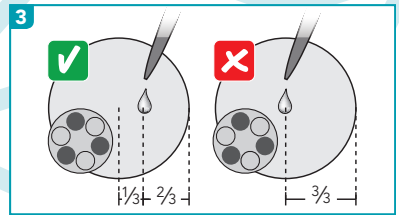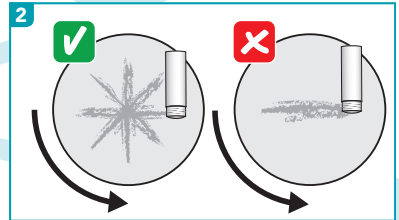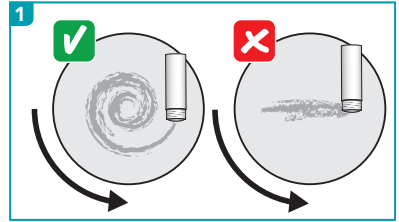

Monokristallin
9 μm - 1 μm
20 ml



Polykristallin
9 μm - 0.25 μm
20 ml

Kontaktieren Sie Ihren ansässigen Akasel Händler um eine bessere Alternative zu Ihrer bisherigen Präparationsmethode zu finden.